



**Ministry of Higher Education
and Scientific Research**





**Al-Mustaqbal University
College of Science**

Department of Medical biotechnology

Biochemistry / Laboratory

Qualitative Analysis of lipids

By:- M.S.c Saja Jawad Abaid

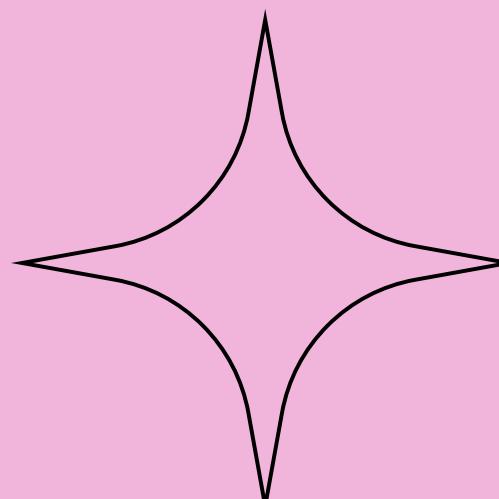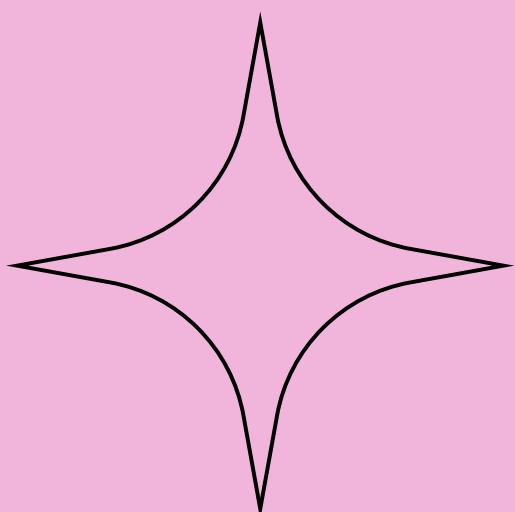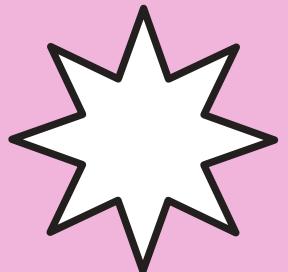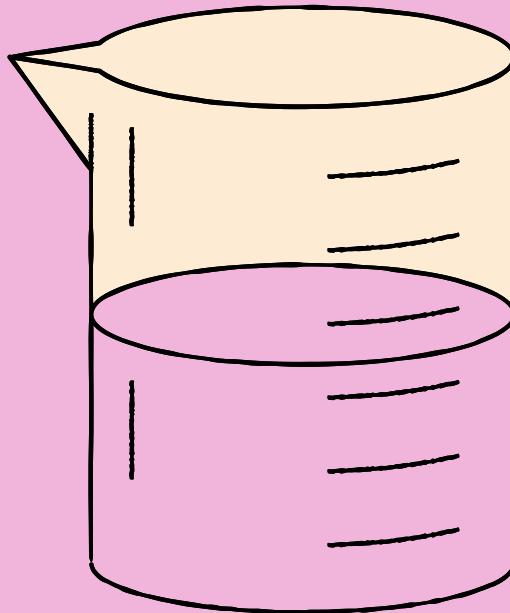




◆ **Definition of Lipids**

Lipids are a group of organic compounds that are insoluble in water but soluble in organic solvents (e.g., ether, chloroform, benzene).

They include fats, oils, waxes, phospholipids, and sterols.

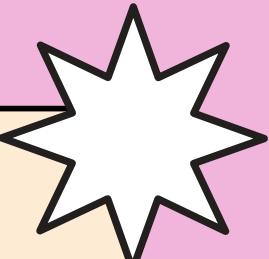

IMPORTANCE OF QUALITATIVE ANALYSIS

- To differentiate between saturated and unsaturated lipids.
- To identify the presence of glycerides and fatty acids.
- To understand lipid reactivity in biological and chemical systems.




Qualitative Tests for Lipids

Test 1: Solubility Test

:Principle

Lipids are nonpolar compounds; hence they are insoluble in water but soluble in organic solvents


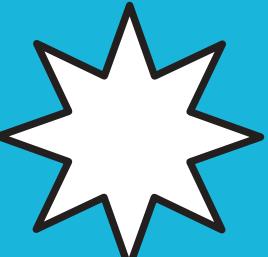


procedure

1 Add 1 mL of the lipid sample into a test tube

2 Add 1 mL of water and shake gently


3 Repeat the test using organic solvents such as ether, chloroform, and ethanol

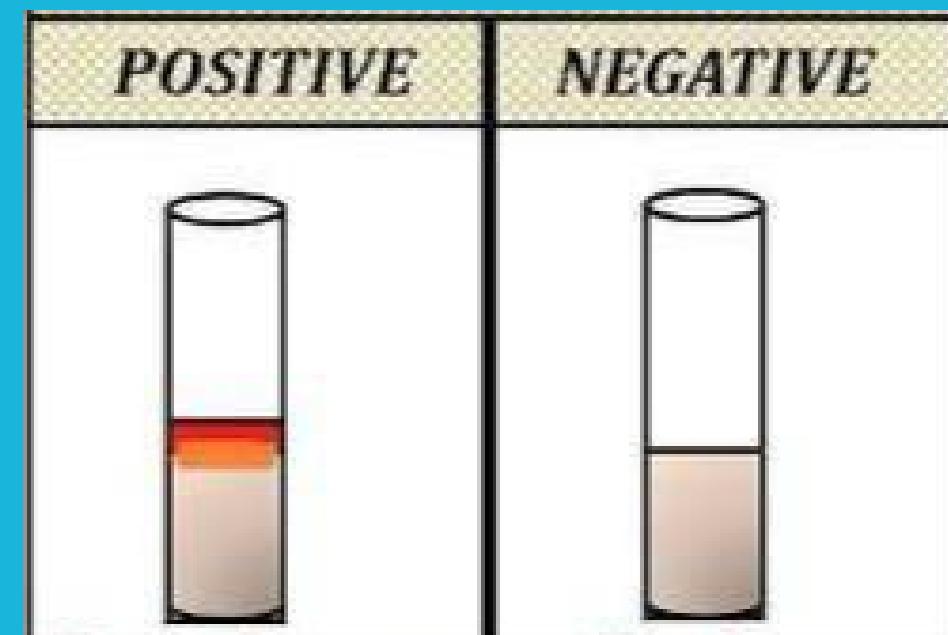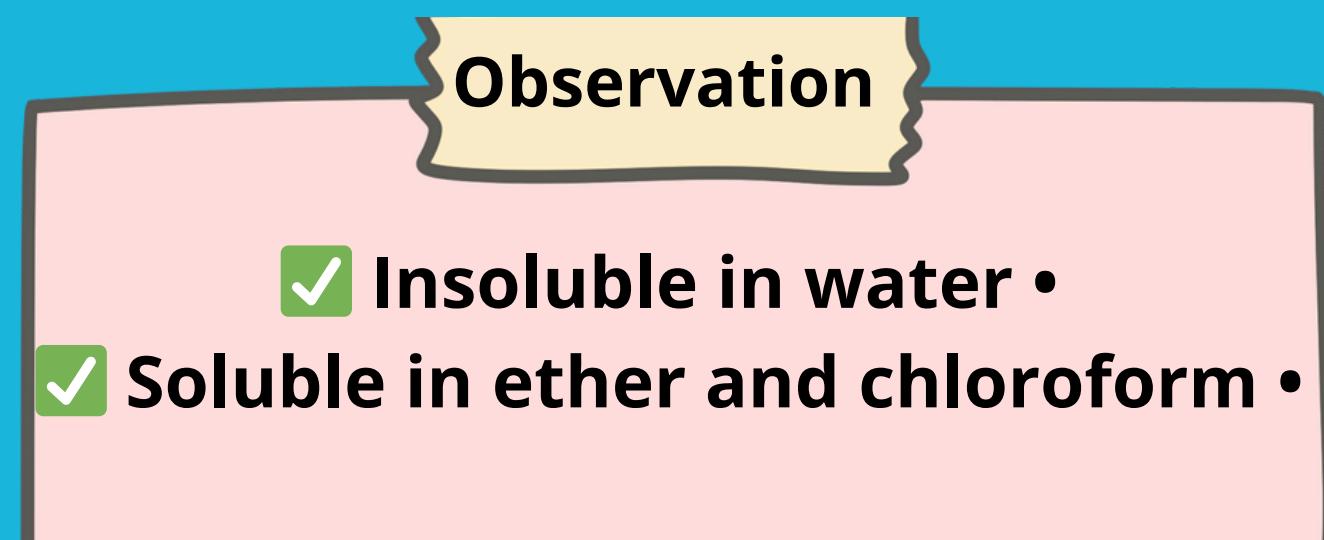

Test 2: Unsaturation Test (Bromine Test)

:Principle

Unsaturated fatty acids contain double bonds that react with bromine, causing the disappearance of bromine's brown color

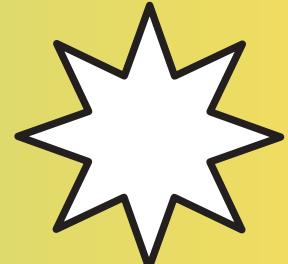

Add a few drops of bromine solution in chloroform to the lipid sample


.Observe the color change

Observation:
Disappearance of • brown color → • Unsaturated lipid
No change in color → • Saturated lipid

Test 3: Acrolein Test (for Glycerides)

:Principle

When glycerides are heated with potassium bisulfate (KHSO_4), acrolein is produced, which has a pungent odor

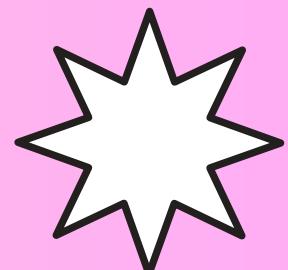

Mix a small amount of lipid sample with KHSO_4 (solid)

Heat gently over a flame





Observation:

- Pungent odor (similar to burnt fat) → Presence of glycerides.**

Test 4: Saponification Test

:Principle

.Alkaline hydrolysis of fats produces soap (fatty acid salt) and glycerol

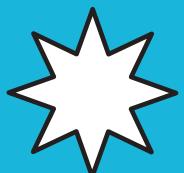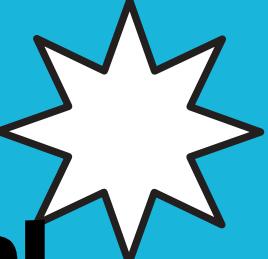

procedure

Add 2 mL of lipid sample to 2 mL
.of alcoholic NaOH solution

Add cold water and
.shake


Observation

- Formation of foam →
Positive test for saponifiable
lipids.

THANK YOU!



”لَا حِيَاةٌ لِمَنْ يَظْلِمُ وَاقْفًا عَلَى
الضَّفَافِ، خَائِفًا مِنَ الْأَمْوَاجِ
وَالْأَعْصِيرِ، الْحِيَاةُ لِمَنْ
يَتَحَركُ، يُقْدِمُ، يُقْبِلُ، يَخُوضُ،
يَتَعَثِّرُ، يَنْهَضُ، يَصْبِرُ، حَتَّى
يَظْفَرُ أَخْيَرًا.”

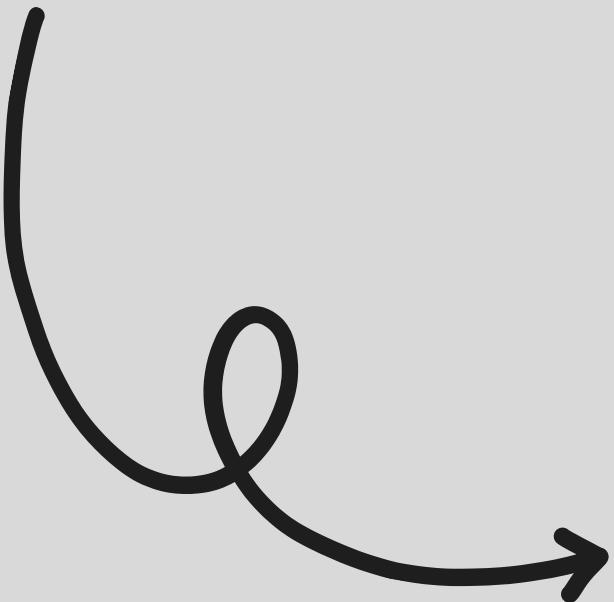

ELEMENTS

